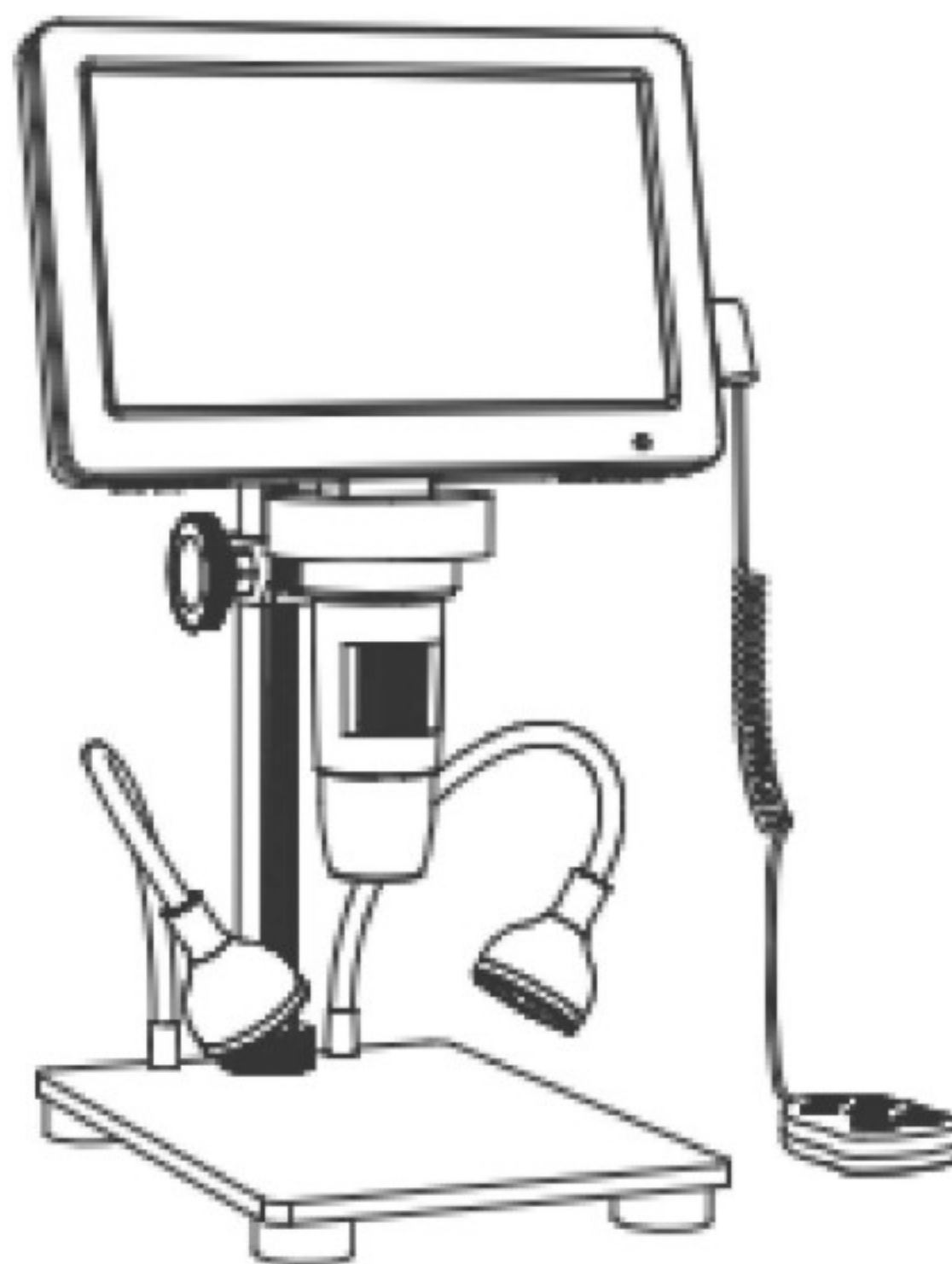
7 Inch Digital Microscope

Model: DM9

Manual


Spec

Power Supply: DC 5V from adapter(need buy it separately), DC 3.7V from the built-in battery	Built-in battery:1800mAh
Charging Time :about 150min	Digital Magnification: 50-1200X
Photo Resolution: 16M, 12M,10M, 8M, 5M, 3M, 2MHD,1.3M	Photo Format: JPG
Video Resolution: 1080FHD,1080P,720P	Video Format: AVI
Light Source: 8 led (brightness adjustable)	Focusing Range: 10~40mm, Manual Focus
PC Compatible With: Windows 7 and above, Mac OSx 12 and above	PC Resolution And Image Transfer Rate: 1280*720/26fps
Support Languages: 13	Operating Temperature: -20°C ~+60°C
Working Humidity: 30%~85%Rh	Power Consumption: 1.95W

INSTALLATION

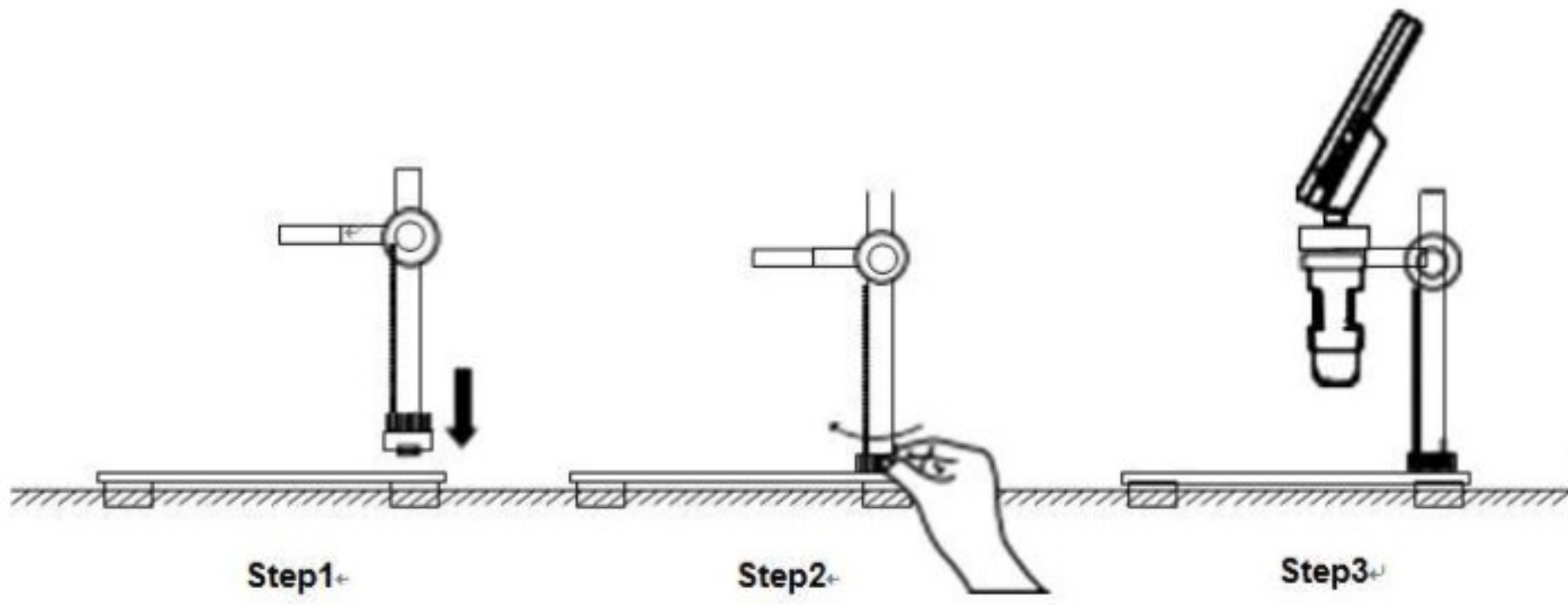1. Metal Stand & Screen Installation

2. Lens Hood Installation

a. Remove the lens cap

b. put on the lens cap

c. Rotate to tighten the lens cap

3. How to connect wired control and filling light cable

PARTS AND FUNCTIONS

Functions introduction

Up/Down Button: it is the selection key.

OK Button:

1. In photo mode, short press this key to take a picture and save it automatically.
2. In video mode, short press it to start/end recording.
3. In playback mode, press it to play the recording.
4. When entering the menu interface, short press this key to confirm the selection.

Photo Button: when in the photo mode, short press it to take a picture.

Menu Button:

1. Short Press this key to enter the menu, and short press this key again to exit the menu and return back to the previous interface.
4. Long press this key to switch between Photo/Video/Playback mode.

Indicator Light: The led is blue after power on; the led will turn from red to green when fully charged.

Power Button: Long press this key for 3 seconds to start/shut down the device.

Type-C Pc Interface: Supports DC 5V recharging and USB data transmission.

(Note: When the USB data cable is removed when link with pc, the device will automatically shut down for protection, which is a normal phenomenon. If you need to continue using the device, you can restart the device.)

TF Slot: Support up to 32G memory card (Note: Please choose **C10 high-speed card**)

Key Slot: wired control and two filling lights cable jacket

Main Led Adjuster: Rotate this wheel to regulate the brightness of the 8 LED.

Focus Wheel: Rotate this wheel to focus on the object.

Reset Button: When the device is abnormal, push this button to restart.

Wired Control Introduction

Switch key: In the power-on status, short press this key to switch between Photo/Video/Playback mode.

Recording key: In recording mode, short press this key to start recording and end recording.

Confirm key: In the photo mode, press this key to take photos and save files automatically. In video mode, short press this key to take video and then press the key to stop and save files. When you enter a menu, you can confirm the selection by short pressing this button.

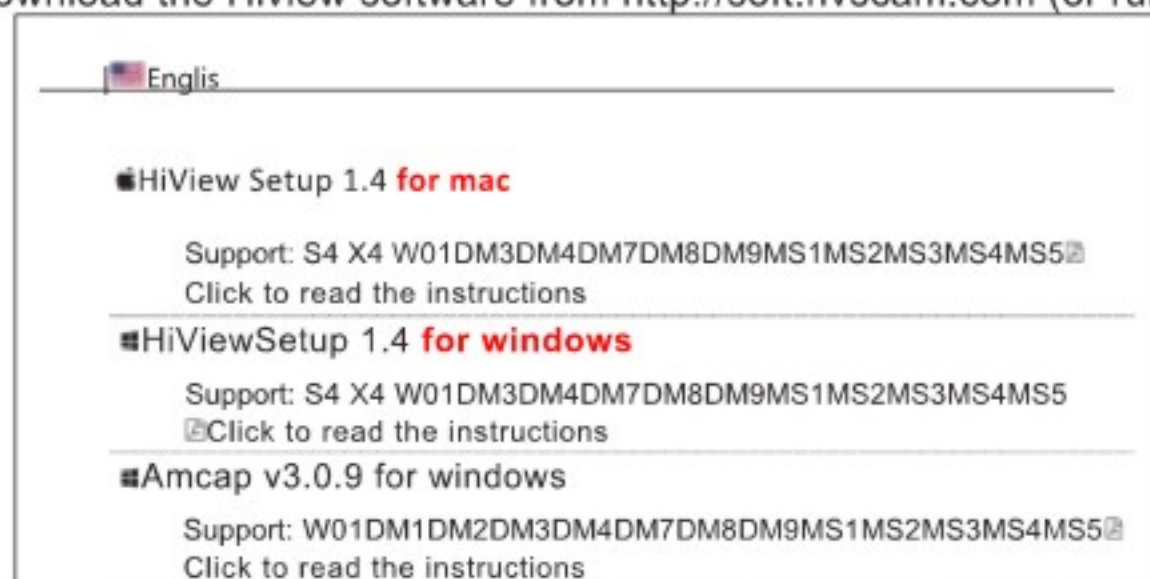
Up/Down key: up /down selection.

Menu key: short press it to enter/exit the menu.

How to Connect With A Pc

Step1

Download the Hiview software from <http://soft.hvscam.com> (or run the software disk to install)


Step 2

Connect the microscope to the PC with a USB cable.

Step 3

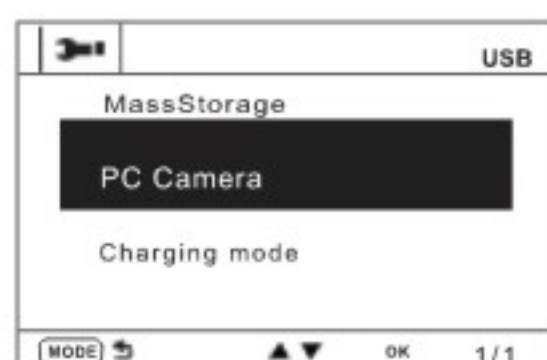Power on DM9, Select "**PC Camera**" on DM9.

Step 4

Run HiView, then you can click Photo or Video Button in Hiview.

Attention

- * To avoid direct sunlight.
- * To avoid heat or fire sources, such as microwave ovens, stoves and other appliances may generate high temperatures.
- * Do not modify the battery, do not immerse it in water
- * Dispose the battery properly, do not throw it into fire may cause explosion

FAQ

Q: Can I see cells by this microscope?

Sorry, this is a digital microscope, not an optical microscope. You are unable to see the cells.

Q : How to recharge it?

There is a USB cable in the package. You can charge the microscope with your phone charger adapter or connect it to your laptop or PC, and the charging time is about 2-5 hours.

Q: When I insert the wire into the key slot, why the microscope show slow battery?

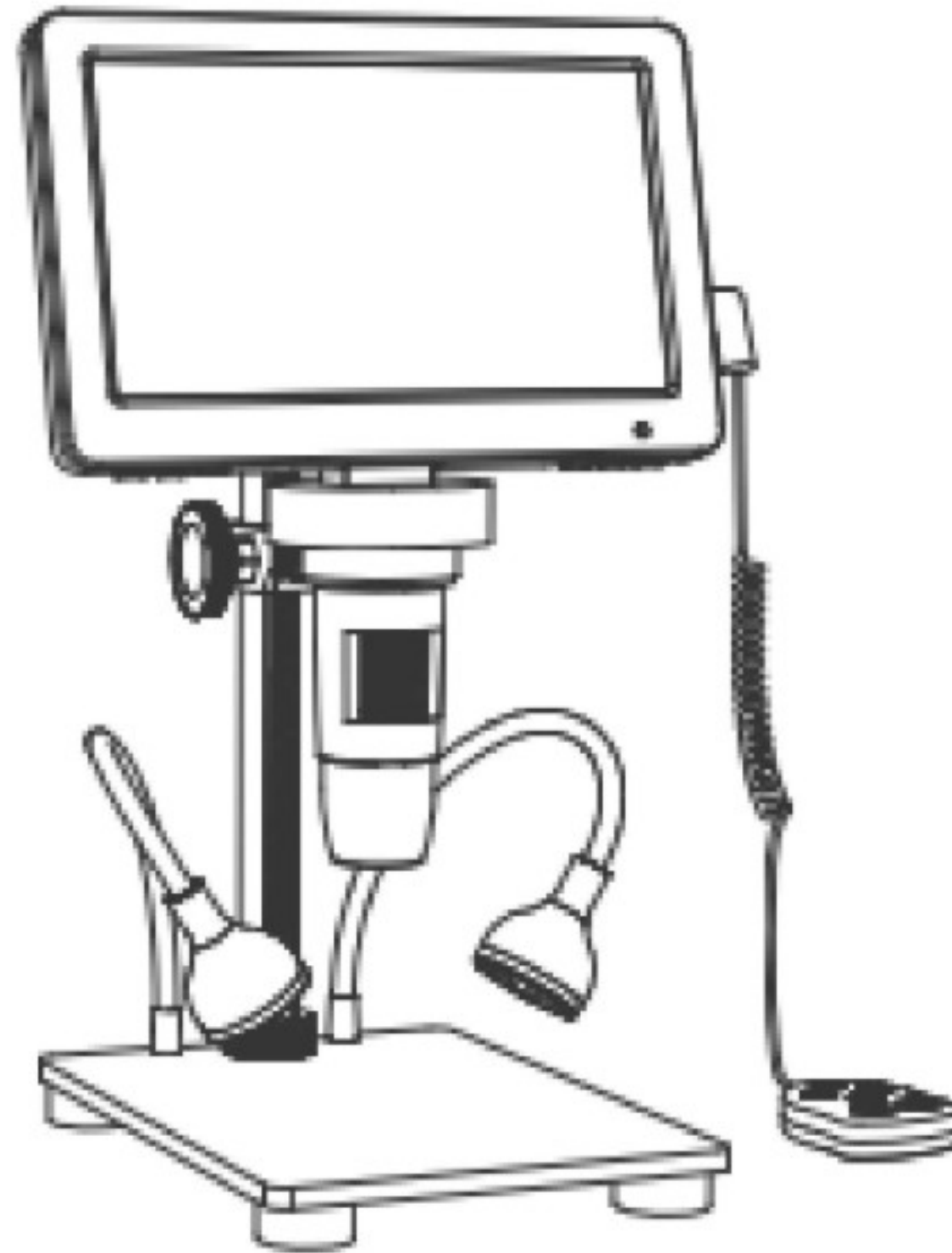
Please insert it into fully position, otherwise it may have the low battery reminding.

Q: Packing List: 1 microscope, 1 metal stand, 1 wired remote control, 1 type cable, 1 lens hood, 1 aid ruler, 1 manual

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